


Visistar™-Integrin Targeted Ultrasound Contrast Agent

Catalog Number: VS-101 1 x 0.3 ml vial


- Ultrasound imaging of angiogenesis
- Targeted to $\alpha_v\beta_3$ integrin
- Ready to administer directly from vial

Applications


Visistar™-Integrin agents can be used for ultrasound molecular imaging of $\alpha_v\beta_3$ integrin expressed on angiogenic vascular endothelium. Agent echogenicity is optimal at ultrasound frequencies in 1-20 MHz range, and ultra-high frequency imaging (20-40 MHz) is also possible¹. Contrast imaging settings such as pulse inversion or CPS must be enabled on the ultrasound scanner for optimal contrast sensitivity. Agents remain acoustically active in vivo for 5-15 minutes, depending on the administered dose and scanner settings. A dose of 30 μ L Visistar-Integrin per 25 g mouse is generally sufficient for tumor imaging; however, dosage should be optimized for each application. Please contact Targeson technical support for protocol assistance.

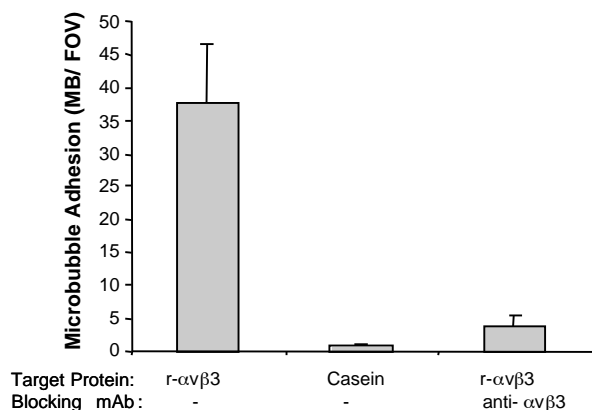

The figure demonstrates adhesion of Visistar-Integrin agents to recombinant mouse $\alpha_v\beta_3$ integrin in a parallel plate flow chamber adhesion assay². The Visistar-Integrin agents exhibit little non-specific adhesion to casein-blocked control surfaces. Incubation of $\alpha_v\beta_3$ flow chambers with an anti- $\alpha_v\beta_3$ blocking antibody abolishes of >95% of Visistar Integrin adhesion.

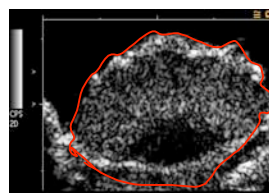

B-mode image

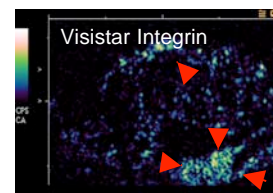

Contrast Image (5 minutes)

Visistar-Integrin in a mouse model of tumor angiogenesis. Human prostate cancer cells (PC-3) implanted subcutaneously on mouse nape, and imaged at day 7. Tumor is outlined in red in the B-mode scan (left). Regions of Visistar-Integrin accumulation are highlighted with red arrows on contrast scan (right). Imaged at 8 MHz using CPS on a Sequoia (Siemens) ultrasound scanner.

Preparation

Visistar-Integrin agents are microbubbles composed of a perfluorocarbon gas core encapsulated by a lipid shell. The agents are further stabilized by a layer of poly(ethylene glycol). The outer shell is derivatized with a peptide that selectively binds endothelial $\alpha_v\beta_3$ integrin. Agents are suspended in aqueous saline at a concentration of approximately 1×10^9 particles per mL, and are packaged in glass vials with a perfluorocarbon gas headspace. The agents have a median diameter of approximately 2.5 μ m.

Storage

Store undiluted at 4°C in sealed vials until ready for application. Concentration is stable in sealed vials for 3 months after septum is punctured. Concentration may decrease upon prolonged storage or vial opening. Do not freeze.

References

1. Rychak JJ, J Graba, AM Cheung, BS Mystry, JR Lindner, RS Kerbel, FS Foster. 2007. Mol Imaging 6(5): 289-96
2. Rychak JJ, B Li, ST Acton, A Leppanen, RD Cummings, K Ley, AL Klibanov. 2006. Mol. Pharm 3(5): 516-24.

FOR RESEARCH USE ONLY. NOT FOR USE IN HUMANS. Targeson makes no representations and extends no warranties of any kind, and expressly disclaims all implied warranties of merchantability and of fitness for a particular purpose or use. Targeson disclaims all warranties of noninfringement with respect to any third party rights and title, including patent rights, in the materials. Use of this product is governed by the Targeson terms and conditions agreement, available at www.targeson.com. Visistar™ is a pending trademark of Targeson Inc.



Targeson, Inc
7590 Fay Ave
Suite 404
La Jolla, CA 92007
Tel: (877) 290-4043
Fax: (877) 401-7805
products@targeson.com
www.targeson.com